
Hygiene Monitoring System

Instructions for use

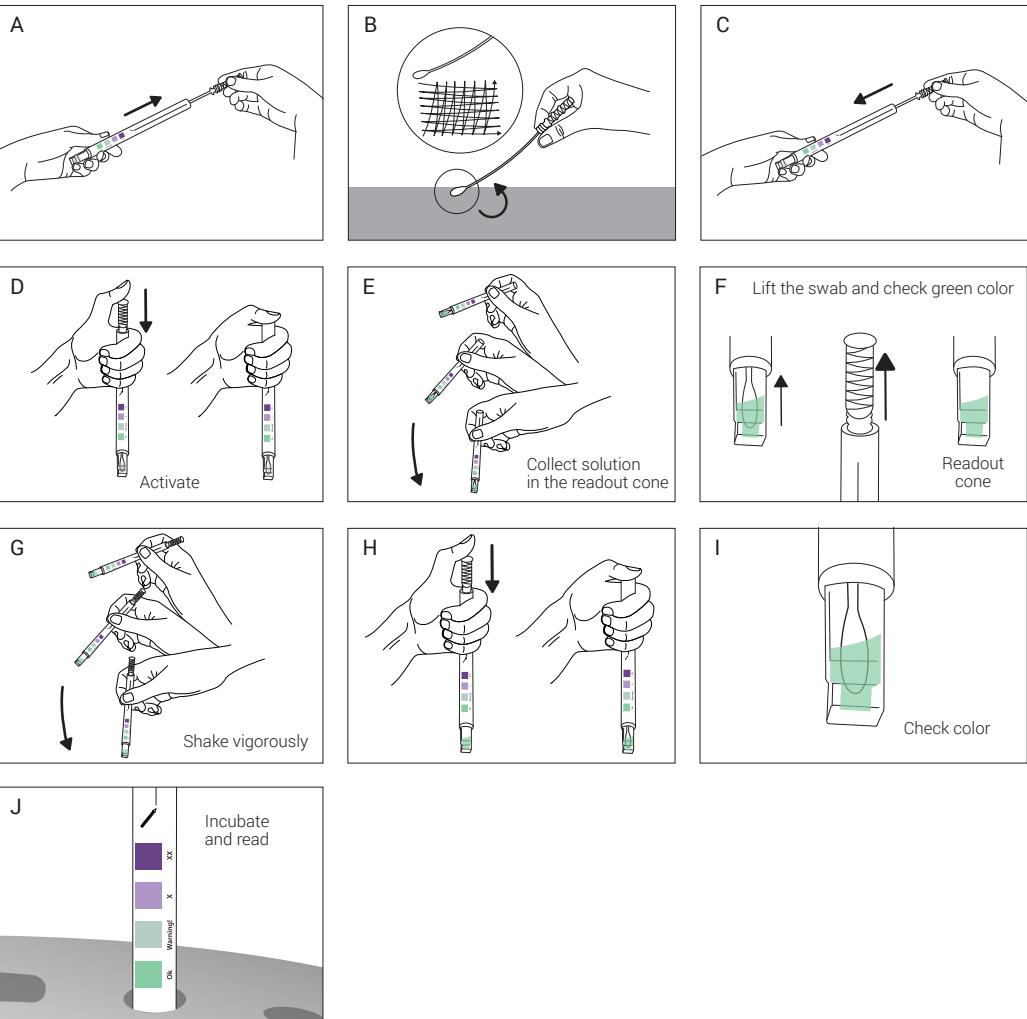PRO1

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PRO 1

Pen system for surface protein and allergen detection

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TR

Product description

Poor cleaning can lead to bacterial growth and furthermore food residues might contain allergens and may represent potential contaminants for allergen-free products. Monitoring the effectiveness of cleaning helps to reduce the risk of a food safety incident.

Protein or Allergen pen system PRO1 has been designed to detect proteins, allergens derived from foods such as milk, egg, gluten, soy and peanut, and to reduce agents on surfaces after the cleaning process and in hard-to-reach areas, in real time. The system has a high absorption swab, that allows collection of samples from different surfaces with the same efficacy.

A visual readout of a color change indicates the presence of detectable levels of protein. PRO1 has high sensitivity and can detect up to 3 µg of protein on surfaces and solutions.

Characteristics

PRO1 provides all the necessary equipment required for HTM2030/HTM01-05 (UK)/ISO15883-1 Residual Protein Test.

Indications for use

Use for testing any surface after the cleaning process and for hard-to-reach areas in complex instruments such as serrated edges or box joints.

Precautions

Do not open until use. Do not reuse. Do not touch the swab with your hands. Wear gloves, surgical mask and other hygiene measures. Do not freeze.

High concentrations of alkaline sanitizers may cause false negative results in some cases. Hydrogen peroxide-based disinfectants may cause a false positive color change reaction.

Although PRO1 has high sensitivity, it is not a microbiological monitoring device, and surfaces may not be considered sterile based on the results obtained.

Instructions for use

1. Remove the swab from the device (fig. A). For dry surfaces, carefully apply 3 drops of moisturizer on the swab or surface to be tested.

2. On the surface to take the sample, swab in one direction and then perpendicularly (fig. B). Surface size should be approximately 10 cm². Apply pressure on the swab and rotate it while you collect the sample.

3. Place the swab back into the device (fig. 1. C) and activate it by pressing it down firmly (fig. D). Shake vigorously downwards for at least 10 seconds (fig. E).

4. In order to correctly mix both solutions contained within the device, it is highly recommended to lift the swab until it is no longer seen within the readout cone (fig. F) and shake downwards vigorously once again for at least 10 seconds (fig. G). The solution will turn to green. Then, press the swab down completely (fig. H) and check color (fig. I).

Caution: Not shaking the device properly so that the liquid of the inner compartment goes down completely, may bring out invalid results.

If you wish to read several swabs simultaneously, once the sample was taken, PRO1 may remain without being activated for a maximum of 4 hours.

5. Incubate the PRO1 during 15 minutes at 55-60 °C or 120 minutes at 37 °C (fig. J).

IMPORTANT: Results analyzed beyond these incubation periods are not valid.

6. Interpret results visually by using the table printed on the label of the device.

By comparing the color of the test against the table an estimation of the surface cleanliness can be made:

OK - Clean, no further action required.

Warning! - Re-clean.

x/xx - Re-clean and re-test.

For a quantitative analysis, you may use Bionova® IC10/20FR Auto Reader. Read the instructions for use of this device previously.

NOTE: For performing comparable cleaning tests, you must standardize surface sampling. In order to achieve this, it is advisable to take the sample always at the same point, with an invariable area of 10 cm², and slide the swab 10 times in one direction and 10 times perpendicularly.

Storage

PRO1 system should be stored at temperatures between 2-25 °C in dark places. Shelf-life under these conditions: 18 months.

Alternatively, if swabs are transported or stored at temperatures not higher than 35 °C, they remain stable for 14 days.

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